



Lupine ELISA Kit

Enzyme Immunoassay for the quantitative determination of Lupine in food

Catalog number: ARG80804

For research use only. Not for use in diagnostic procedures.

TABLE OF CONTENTS

SECTION	Page
INTRODUCTION	3
PRINCIPLE OF THE ASSAY	4
MATERIALS PROVIDED & STORAGE INFORMATION.....	4
MATERIALS REQUIRED BUT NOT PROVIDED.....	5
TECHNICAL HINTS AND PRECAUTIONS	5
SAMPLE COLLECTION & STORAGE INFORMATION	5
REAGENT PREPARATION	6
ASSAY PROCEDURE	7
CALCULATION OF RESULTS.....	7
EXAMPLE OF TYPICAL STANDARD CURVE.....	8
QUALITY ASSURANCE	9

MANUFACTURED BY:

Arigo Biolaboratories Corporation

Address: No. 22, Ln. 227, Gongyuan Rd., Hsinchu City 300, Taiwan

Phone: +886 (3) 562 1738

Fax: +886 (3) 561 3008

Email: info@arigobio.com

INTRODUCTION

Lupin belongs to the legume family. With about 35-45% the fraction of proteins in lupin seed is very high. Some of these proteins, predominantly those in the α -, β - and γ -Conglutin fractions are known for being allergenic, due to cross-reaction to allergens of soy and peanut. Because of its convenient nutritional properties the use of lupin flour as a food ingredient is highly increasing. Lupin is often used in dietary products, especially for milk allergic people, vegetarians and patients with celiac disease. In food production three (sweet) lupin species are commonly used: *Lupinus albus*, *Lupinus angustifolius* and *Lupinus luteus*.

For lupin allergic persons hidden lupin allergens in food are a critical problem. Already very low amounts of lupin can cause allergic reactions, which may lead to anaphylactic shock in severe cases. Because of this, lupin allergic persons must strictly avoid the consumption of lupin containing food. Cross-contamination, mostly in consequence of the production process, is often noticed. This explains why in many cases the existence of lupin residues in food cannot be excluded. For this reason sensitive detection systems for lupin residues in food-stuffs are required.

The Lupin ELISA represents a highly sensitive detection system for lupin and is particularly capable of the quantification of residues in sausage, bakery products, potato products, ketchup and juices. Due to high cross-reactivity the test is suitable for the detection of all commonly used lupin species.

Lupine ELISA Kit ARG80804

PRINCIPLE OF THE ASSAY

This assay employs the sandwich quantitative enzyme immunoassay technique. An antibody directed against Lupine is bound on the surface of a microtiter plate. Lupine containing samples or standards are given in-to the wells of the microtiter plate. After 20 minutes incubation at room temperature, the wells are washed with diluted washing solution to remove unbound material. A peroxidase conjugated second antibody directed against Lupine proteins is given into the wells and after 20 minutes of incubation the plate is washed again. A substrate solution is added and incubated for 20 minutes, resulting in the development of a blue color. The color development is inhibited by the addition of a stop solution, and the color turns yellow. The yellow color is measured at 450 nm. The concentration of Lupine is directly proportional to the color intensity of the test sample.

MATERIALS PROVIDED & STORAGE INFORMATION

Store the unopened kit at 2-8 °C. Use the kit before expiration date.

Component	Quantity	Storage information
Antibody-coated microplate	12 strips x 8-well	4°C
HRP-antibody Conjugate	15 ml (ready to use)	4°C
Standards (0,2, 5, 15, 30 ppm)	5 X 1 ml (ready to use)	4°C
10x Extraction and sample dilution buffer	2 X 120 ml	4°C
10x Wash Buffer	60 ml	4°C
TMB substrate	15 ml	4°C (Protect from light)
STOP solution	15 ml	4°C

MATERIALS REQUIRED BUT NOT PROVIDED

- Microplate reader capable of measuring absorbance at 450nm
- Pipettes and pipette tips
- Deionized or distilled water
- Automated microplate washer (optional)

TECHNICAL HINTS AND PRECAUTIONS

- Wear protective gloves, clothing, eye, and face protection especially while handling blood or body fluid samples.
- Store the kit at 4°C at all times.
- Briefly spin down the HRP-Antibody conjugate before use.
- If crystals are observed in the 10X Wash buffer, Extraction Buffer and Sample diluent buffer, warm to RT (not more than 50°C) until the crystals are completely dissolved.
- Ensure complete reconstitution and dilution of reagents prior to use.
- It is highly recommended that the standards, samples and controls be assayed in duplicates.
- Change pipette tips between the addition of different reagent or samples.
- Samples contain azide cannot be assayed.

SAMPLE COLLECTION & STORAGE INFORMATION

Due to high risk of cross-contamination all applied instruments like applicator, mortar, glass vials etc. have to be cleaned thoroughly before and after each sample. Hazelnut proteins adhere very strongly to different surfaces. In

Lupine ELISA Kit ARG80804

certain cases they can resist a common dishwasher cleaning. To identify possible cross-contamination caused by previous extractions it is strongly recommended to note the sequence of the extractions.

The following sample preparation should be applied for all kinds of samples:

1. To maximize homogeneity and representativeness of the sample drawing, a minimum of 5 g sample should be pulverized finely in a mortar, impact mill etc.
2. 1 g of the homogenized mixture is suspended in 20 mL of pre-diluted extraction buffer. Afterwards the suspension is incubated for 15 min in a preheated water bath at 60°C. To ensure good homogeneity, the samples should be shaken every two minutes.
3. The samples are centrifuged for 10 minutes at 2000 g. If it is not possible to separate the supernatant from the precipitate completely, the suspension should be filtrated if necessary.
4. 100 µL of particle-free solution are applied per well. If the results of a sample are out of the measuring range, further dilution with the pre-diluted extraction and sample dilution buffer is necessary. The additional dilution has to be considered when calculating the concentration.

REAGENT PREPARATION

- **1X Wash buffer:** Dilute 10X wash buffer into distilled water to yield 1X wash buffer.
- **1X Extraction and Sample diluent buffer:** Dilute 10X Extraction and

Sample diluent buffer into distilled water to yield 1X.

ASSAY PROCEDURE

All materials should be equilibrated to room temperature (RT) before use. Standards, samples and controls should be assayed in duplicates.

1. Remove excess microplate strips from the plate frame, return them to the foil pouch containing the desiccant pack, and reseal it.
2. Add 100 μ l of standards and samples in duplicate into wells.
3. Incubate for 20 minutes at RT.
4. Aspirate each well and wash, repeating the process 2 times for a total 3 washes. Wash by filling each well with 1X wash buffer (350 μ l) using a squirt bottle, manifold dispenser, or autowasher. Complete removal of liquid at each is essential to good performance. After the last wash, remove any remaining Wash Buffer by aspirating, decanting or blotting against clean paper towels.
5. Add 100 μ l of HRP-Antibody Conjugate into each well. Incubate for 20 minutes at RT.
6. Aspirate and wash well as step 4.
7. Add 100 μ l of TMB mixture to each well. Incubate for 20 minutes at room temperature in dark.
8. Add 100 μ l of Stop Solution to each well.
9. Read the OD with a microplate reader at 450 nm immediately.

CALCULATION OF RESULTS

1. Calculate the average absorbance values for each set of standards, controls


Lupine ELISA Kit ARG80804

and patient samples.

2. Using linear graph paper, construct a standard curve by plotting the mean absorbance obtained from each standard against its concentration with absorbance value on the vertical (Y) axis and concentration on the horizontal (X) axis.
3. Using the mean absorbance value for each sample determine the corresponding concentration from the standard curve.
4. Automated method: The results in the IFU have been calculated automatically using a 4 PL (4 Parameter Logistics) curve fit. 4 Parameter Logistics is the preferred method. Other data reduction functions may give slightly different results. The diluted samples must be further converted by the appropriate dilution factor according to the sample preparation procedure as described above.

EXAMPLE OF TYPICAL STANDARD CURVE

The following data is for demonstration only and cannot be used in place of data generations at the time of assay.

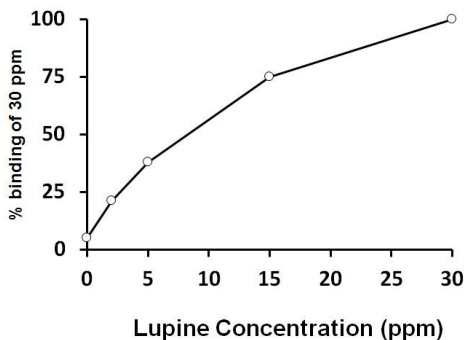

QUALITY ASSURANCE

Sensitivity

The limit of detection (LOD) of the Lupine test is 0.2 ppm for the standard curve.

Validation experiments with common matrices resulted in the following LODs [ppm].

Sausage	0.2
Bread	0.3
Orange juice	0.7
Ketchup	0.1
Croquette	0.2

The limit of quantification (LOQ) of the Lupine test is 2 ppm.

Due to the variety of sample matrices and their influence on the blank, results less than the LOQ should be treated as negative.

Specificity

For the following foods no cross-reactivity could be detected:

Milk	Sunflower seeds	Cocoa
Egg	Pumpkin seeds	Potato
Wheat	Cashew	Leek
Rye	Peanut	Pea
Oats	Hazelnut	Bean
Barley	Almond	Tomato
Rice	Pecan	Peach
Corn	Coconut	Cherry
Buckwheat	Brazil nut	Orange
Sesame	Walnut	Strawberry
Pork	Pistachio	Plum
Beef	Macadamia nut	Locats bean gum
Chicken	Chestnut	Guar gum

Lupine ELISA Kit ARG80804

The following cross reactions were determined:

Chickpea	0.0003%
Lentil	0.0004%
Soy flour, unroasted	0.07%
Soy flour, roasted	0.0009%

Intra-assay and Inter-assay precision

The CV value of intra-assay precision was 6-9% and the CV value of inter-assay precision was 5-9%

Recovery

Sausage	99%
Bread	113%
Orange juice	104%
Ketchup	98%
Croquette	111%