



HSP70 (High sensitive) ELISA Kit

Enzyme Immunoassay for the quantitative determination of HSP70 in cell/tissue lysates, plasma and serum.

Catalog number: ARG81149

For research use only. Not for use in diagnostic procedures.

TABLE OF CONTENTS

SECTION	Page
INTRODUCTION	3
PRINCIPLE OF THE ASSAY	3
MATERIALS PROVIDED & STORAGE INFORMATION.....	3
MATERIALS REQUIRED BUT NOT PROVIDED	5
TECHNICAL HINTS AND PRECAUTIONS	5
SAMPLE COLLECTION & STORAGE INFORMATION	7
REAGENT PREPARATION.....	7
ASSAY PROCEDURE	8
CALCULATION OF RESULTS	10
EXAMPLE OF TYPICAL STANDARD CURVE	11
QUALITY ASSURANCE.....	11

MANUFACTURED BY:

Arigo Biolaboratories Corporation

Address: No. 22, Ln. 227, Gongyuan Rd., Hsinchu City 300, Taiwan

Phone: +886 (3) 562 1738

Fax: +886 (3) 561 3008

Email: info@arigobio.com

INTRODUCTION

HSP70 is a 70kDa heat shock protein which is a member of the heat shock protein 70 family. In conjunction with other heat shock proteins, this protein stabilizes existing proteins against aggregation and mediates the folding of newly translated proteins in the cytosol and in organelles. It is also involved in the ubiquitin-proteasome pathway through interaction with the AU-rich element RNA-binding protein 1. The gene is located in the major histocompatibility complex class III region, in a cluster with two closely related genes which encode similar proteins. [provided by RefSeq, Jul 2008]

HSP70 is in cooperation with other chaperones, Hsp70s stabilize preexistent proteins against aggregation and mediate the folding of newly translated polypeptides in the cytosol as well as within organelles. These chaperones participate in all these processes through their ability to recognize nonnative conformations of other proteins. They bind extended peptide segments with a net hydrophobic character exposed by polypeptides during translation and membrane translocation, or following stress-induced damage. In case of rotavirus A infection, serves as a post-attachment receptor for the virus to facilitate entry into the cell. Essential for STUB1-mediated ubiquitination and degradation of FOXP3 in regulatory T-cells (Treg) during inflammation. [UniProt]

PRINCIPLE OF THE ASSAY

This assay employs the quantitative sandwich enzyme immunoassay technique. An antibody specific for HSP70 has been pre-coated onto a microtiter plate. Standards or samples are pipetted into the wells and any HSP70 present is

HSP70 (High sensitive) ELISA Kit ARG81149

bound by the immobilized antibody. After washing away any unbound substances, a biotin-conjugated antibody specific for HSP70 is added to each well and incubate. Following a washing to remove unbound substances, streptavidin conjugated to Horseradish Peroxidase (HRP) is added to each microplate well and incubated. After washing away any unbound antibody-enzyme reagent, a substrate solution (TMB) is added to the wells and color develops in proportion to the amount of HSP70 bound in the initial step. The color development is stopped by the addition of acid and the intensity of the color is measured at a wavelength of 450nm \pm 2nm. The concentration of HSP70 in the sample is then determined by comparing the O.D of samples to the standard curve.

MATERIALS PROVIDED & STORAGE INFORMATION

Store the unopened kit at 4°C. Use the kit before expiration date.

Component	Quantity	Storage information
Antibody-coated microplate	96 wells	4°C
Standard (Lyophilized)	2 vials	4°C
Standard/Sample diluent	50 ml (ready to use)	4°C
100X Antibody conjugate concentrate	150 μ l	4°C
Antibody diluent buffer	13 ml (ready to use)	4°C
500X HRP-Streptavidin concentrate	50 μ l	4°C
HRP-Streptavidin diluent buffer	13 ml (ready to use)	4°C
10X Wash buffer	100 ml	4°C

HSP70 (High sensitive) ELISA Kit ARG81149

TMB substrate	13 ml (ready to use)	4°C (Protect from light)
STOP solution	13 ml (ready to use)	4°C

MATERIALS REQUIRED BUT NOT PROVIDED

- Microplate reader capable of measuring absorbance at 450 nm
- Pipettes and pipette tips
- Deionized or distilled water
- Polypropylene or polyethylene tubes to prepare samples – do not use polystyrene, polycarbonate or glass tubes.
- Adhesive plate sealers
- Automated microplate washer (optional)
- Tissue/Cell protein Extraction Reagent (Not Supplied, optional):
 - a. Extraction Reagents must contain $\leq 0.1\%$ SDS eg. 50mM Tris pH7.2, 150 mM NaCl, 1% Triton X-100.

Note: Alternative extraction reagents may be used, but reagents containing more than 0.1% SDS will affect results.

- b. Add protease inhibitors to the Extraction Reagent before use. Examples of an appropriate protease inhibitor cocktail includes 0.1 mM PMSF, 1 $\mu\text{g}/\text{mL}$ leupeptin, 1 $\mu\text{g}/\text{mL}$ aprotinin, and 1 $\mu\text{g}/\text{mL}$ pepstatin. Alternatively, a commercially available protease cocktail, obtainable from a variety of scientific reagent vendors, may also be used.

TECHNICAL HINTS AND PRECAUTIONS

- Wear protective gloves, clothing, eye, and face protection especially while handling blood or body fluid samples.
- Store the kit at 4°C at all times.
- If crystals are observed in the 10X Wash buffer, warm to RT (not more than 50°C) until the crystals are completely dissolved.
- Ensure complete reconstitution and dilution of reagents prior to use.
- All materials should be equilibrated to room temperature (RT, 22-25°C) 20 min before use.
- All reagents should be mixed by gentle inversion or swirling prior to use. Do not induce foaming.
- Before using the kit, spin tubes and bring down all components to the bottom of tubes.
- Mix the contents of the microplate wells thoroughly by microplate shaker for few seconds or gently tap the plate to ensure good test results. Please mix carefully to avoid well-to-well contamination. Do not reuse microwells.
- The TMB substrate should be colorless and transparent before using. Take care not to contaminate the TMB Substrate. Do not expose TMB Substrate solution to glass, foil, or metal. If the solution is blue before use, DO NOT USE IT.
- Use reservoirs only for single reagents. This especially applies to the substrate reservoirs. Using a reservoir for dispensing a substrate solution that had previously been used for the conjugate solution may turn

HSP70 (High sensitive) ELISA Kit ARG81149

solution colored. Do not pour reagents back into vials as reagent contamination may occur.

- Do not let wells dry during assay; add reagents immediately after completing the rinsing steps.
- Avoid using reagents from different batches.
- It is highly recommended that the standards, samples and controls be assayed in duplicates.

Change pipette tips between the addition of different reagent or samples.

SAMPLE COLLECTION & STORAGE INFORMATION

The sample collection and storage conditions listed below are intended as general guidelines. Sample stability has not been evaluated.

Serum- Use a serum separator tube (SST) and allow samples to clot for 30 minutes before centrifugation for 10 minutes at 2400 x g. Collect serum and assay immediately or aliquot and store samples at $\leq -20^{\circ}\text{C}$. Avoid repeated freeze-thaw cycles.

Plasma - Collect plasma using EDTA, heparin or Sodium Citrate as an anticoagulant. Centrifuge for 10 minutes at 2400 x g within 30 minutes of collection. Assay immediately or aliquot and store samples at $\leq -20^{\circ}\text{C}$. Avoid repeated freeze-thaw cycles.

Cell Lysate Preparation-

1. Prepare and treat cells as desired.
2. a. For adherent cells- Remove media and rinse cells with ice-cold PBS. Harvest cells with trypsin-EDTA or by using a cell scraper. Centrifuge at 500 x g for 5 minutes.

HSP70 (High sensitive) ELISA Kit ARG81149

- b. For suspension cells- Harvest by centrifugation at 500 x g for 5 minutes.
3. Wash cells by re-suspending the cell pellet in ice-cold PBS. Pellet cells by centrifugation at 500 x g for 5 minutes. Repeat wash for a total of three washes with ice-cold PBS.
4. Use a pipette to carefully remove and discard the supernatant, leaving the cell pellet as dry as possible. The cell pellet may be frozen at -70°C and lysed at a later time, if desired.
5. Calculate the amount of Extraction Reagent required (Not Supplied).
For every 1×10^6 to 1×10^7 cells, use 1 mL of 1X Extraction Reagent.
Note: Alternative extraction reagents may be used, but reagents containing more than 0.1% SDS will affect results
6. Add protease inhibitors to the Extraction Reagent before use.
7. Add appropriate amount of ice-cold Extraction Reagent including protease inhibitors to the cell pellet.
Note: If excess buffer is used for the number of cells lysed, the protein concentration will be low.
8. Pipet up and down to break up the cell pellet until the cell suspension is homogeneous and no clumps are visible.
9. Incubate on ice for 30 minutes with occasional mixing or sonication.
Note: To increase protein yields and decrease sample viscosity, aspirate the cell pellet 5-10 times through a 21 ½ gauge needle or sonicate the cell pellet for 30 seconds with 50% pulse during the incubation.
10. Transfer the mixture to a fresh micro centrifuge tube and centrifuge at ~21,000 x g for 10 minutes at 4°C.

HSP70 (High sensitive) ELISA Kit ARG81149

11. Transfer the supernatant (cell lysate) to a fresh tube for analysis. Avoid disturbing the cell pellet. Discard the cell pellet once the supernatant is harvested. The cell lysate is now ready for analysis in the assay.
12. Alternatively, store the cell lysate in single-use aliquots at -70°C . It is recommended that a protein determination assay be performed and the extracts aliquoted into convenient amounts prior to storing at -70°C to avoid multiple freeze-thaw cycles.

Tissue Extract Preparation-

1. Harvest tissue to be analyzed. Tissues may be flash frozen, stored at -70°C and prepared at a later time, if desired.
2. Calculate the amount of Extraction Reagent required (Not Supplied). For each 0.5 cm^3 piece of tissue, use 1 mL of Extraction Reagent.

Note: Alternative extraction reagents may be used, but reagents containing more than 0.1% SDS will affect results.

3. Add protease inhibitors to the Extraction Reagent.
4. Place the tissue in a mortar and add sufficient volume of liquid nitrogen to cover the tissue.
5. Allow the liquid nitrogen to evaporate. The tissue should be thoroughly frozen.
6. Grind the frozen tissue to a powder with a pestle.
7. Add appropriate amount of ice-cold Extraction Reagent including protease inhibitors to the processed tissue. Continue to homogenize the tissue with the pestle until the tissue suspension is homogeneous.
9. Transfer the extract to a fresh micro centrifuge tube and centrifuge at $\sim 21,000 \times g$ for 10 minutes at 4°C .

HSP70 (High sensitive) ELISA Kit ARG81149

10. Transfer the supernatant (tissue extract) to a fresh tube for analysis. Avoid disturbing the cell pellet. Discard the cell pellet once the supernatant is harvested. The tissue extract is now ready for analysis in the assay.
11. Alternatively, store the tissue extract in single-use aliquots at -70°C . It is recommended that a protein determination assay be performed and the extracts aliquoted into convenient amounts prior to storing at -70°C to avoid multiple freeze-thaw cycles.

Note:

- a) Do not use haemolytic, icteric or lipaemic specimens.
- b) Samples containing sodium azide should not be used in the assay.
- c) The cell lysate Extraction Reagent is not included with this kit.
- d) If particulate is present in samples, centrifuge prior to analysis.

REAGENT PREPARATION

- **1X Wash buffer:** Dilute 10X Wash buffer into distilled water to yield 1X Wash buffer. (E.g. 100 ml of 10X Wash buffer + 900 ml of distilled water)
The diluted Wash buffer is stable for one month at 4°C. Do not use 1X Wash Buffer if it becomes visibly contaminated during storage.
- **Antibody conjugate working solution:** Prior to use, prepare only the required amount of 1X Antibody conjugate working solution for the number of strips being used. Dilute Antibody conjugate concentrate (100X) into Antibody diluent buffer and mix well to yield 1X Antibody conjugate working solution. (E.g. 6 ml of 1X Antibody conjugate working solution is required for 48 wells used: 60 µl of Antibody conjugate concentrate (100X) + 6 ml of Antibody diluent buffer, mix well)
- **HRP-Streptavidin working solution:** Prior to use, prepare only the required amount of 1X HRP-Streptavidin working solution for the number of strips being used. Dilute HRP-Streptavidin concentrate (500X) into HRP-Streptavidin diluent buffer and mix well to yield 1X HRP-Streptavidin conjugate working solution. (E.g. 6 ml of 1X HRP-Streptavidin working solution is required for 48 wells used: 12 µl of HRP-Streptavidin conjugate concentrate (500X) + 6 ml of HRP-Streptavidin diluent buffer, mix well)
- **Sample Dilution:** Samples must first be diluted prior to testing.
 - Suggested starting dilutions for samples:
 - For native human cell lysates, dilute samples 1:128 in Standard/Sample diluent. For example, dilute 2 µL of sample in 254

HSP70 (High sensitive) ELISA Kit ARG81149

μ L Standard/Sample diluent. Mix well. For non-human cell lysates, dilute samples 1:4 in Standard/Sample diluent.

- For serum and plasma samples, dilute samples 1:4 in Standard/Sample diluent. For example, dilute 60 μ L of sample in 180 μ L Standard/Sample diluent. Mix well.

Note: If values for samples are not within the range of the standard curve, optimal sample dilutions need to be determined.

- Prepare at least 240 μ L of sample in Standard/Sample diluent. Mix samples well prior to analysis.
- **Standards:** Reconstitute the standard with 1 ml Standard/Sample diluent to yield a stock concentration of 54 ng/ml. Keep the buffer in the vial for at least 5 min at RT to make sure the standard is dissolved completely before making serial dilutions. The standard diluent buffer serves as zero standard (0 ng/ml), and the rest of the standard serial dilution can be diluted as according to the suggested concentration below: 27 ng/ml, 13.5 ng/ml, 6.75 ng/ml, 3.38 ng/ml, 1.69 ng/ml, 0.84 ng/ml, 0.42 ng/ml.

Dilute HSP70 standard as according to the table below:

Standard	HSP70 Conc. (ng/ml)	μ l of Assay Buffer	μ l of standard
S7	27	500	500 (54 ng/ml stock)
S6	13.5	250	250 (S7)
S5	6.75	250	250 (S6)
S4	3.38	250	250 (S5)
S3	1.69	250	250 (S4)
S2	0.84	250	250 (S3)
S1	0.42	250	250 (S2)
S0	0	250	0

ASSAY PROCEDURE

All materials should be equilibrated to room temperature (RT) before use. Standards, samples and controls should be assayed in duplicates.

1. Remove excess microtiter strips from the plate frame, return them to the foil pouch containing the desiccant pack, and reseal it.
2. Add 100 μ l of the Standards and diluted samples into the appropriate wells. Gently tap the plate to mix well. Cover the plate and incubate for 2 h at 37°C.
3. Aspirate each well and wash, repeating the process 3 times for a total 4 washes. Wash by filling each well with 1 \times Wash Buffer (300 μ l) using a squirt bottle, manifold dispenser, or autowasher. Complete removal of liquid at each is essential to good performance. After the last wash, remove any remaining Wash Buffer by aspirating, decanting or blotting.
4. Add 100 μ l of the 1X Antibody conjugate working solution into the appropriate wells. Gently tap the plate to mix well. Cover the plate and incubate for 2 h at 37°C.
5. Aspirate each well and wash as step 3.
6. Add 100 μ l of the 1X HRP-Streptavidin working solution into the appropriate wells. Gently tap the plate to mix well. Cover the plate and incubate for 30 min at 37°C.
7. Aspirate each well and wash as step 3.
8. Add 100 μ l of TMB substrate solution into each well. Incubate for 30 mins at RT without shaking. Avoid exposure to light.
9. Add 100 μ l of Stop Solution to each well and shake or tap the plate lightly

HSP70 (High sensitive) ELISA Kit ARG81149

to ensure homogeneous mixing. The color of the solution should change from blue to yellow.

10. Read the OD with a microplate reader at 450nm immediately. It is recommended read the absorbance within 30 min after adding Stop solution.


CALCULATION OF RESULTS

1. Calculate the average absorbance values for each set of standards, controls and patient samples.
2. Using linear graph paper, construct a standard curve by plotting the mean absorbance obtained from each standard against its concentration with absorbance value on the vertical (Y) axis and concentration on the horizontal (X) axis.
3. Using the mean absorbance value for each sample determine the corresponding concentration from the standard curve.
4. Automated method: The results in the IFU have been calculated automatically using a 4 PL (4 Parameter Logistics) curve fit. 4 Parameter Logistics is the preferred method. Other data reduction functions may give slightly different results.
5. If the samples have been diluted, the concentration read from the standard curve must be further converted by the appropriate dilution factor according to the sample preparation procedure as described above.

HSP70 (High sensitive) ELISA Kit ARG81149

EXAMPLE OF TYPICAL STANDARD CURVE

The following data is for demonstration only and cannot be used in place of data generations at the time of assay.

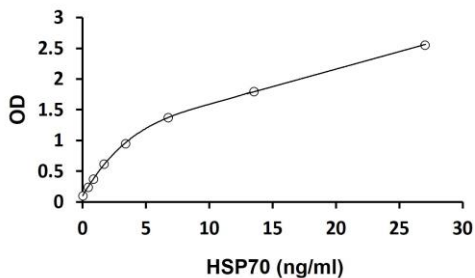

QUALITY ASSURANCE

Sensitivity

The minimum detectable dose (MDD) of HSP70 ranged from 0.42-27 ng/ml. The mean MDD was 0.031 ng/ml.

Specificity

This ELISA kit is specific for Hsp70. This HSP70 ELISA kit has been validated for the detection of Hsp70 from Human, Monkey, Canine, Rat, and Mouse.